


Title:	BCR Bad Reading at Transfer Station			ID:
				0010
Date in:	Response:	Model:	Author:	
2006-08-14	2006-08-14	-	CMa	

Q:

We observe unreliable reading of the BCR at the transfer station. The beam is centered nicely on the barcode. The reading in the stackers is ok.

A:

On clear plates there could be interferences between the BCR and the plate detection sensor of the transfer station.

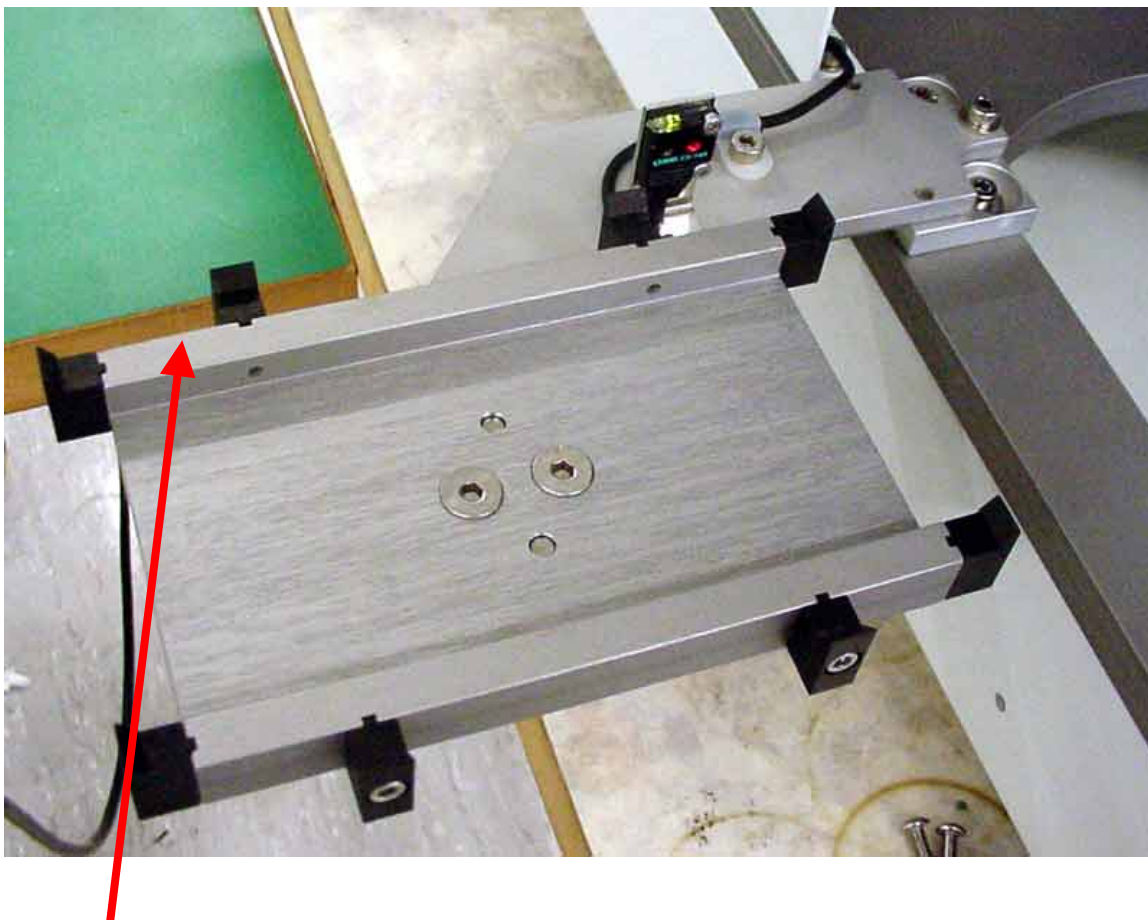

Please have the sensor placed in the location of the Arrow, this is what I did for Astra on the STX 1000 to eliminate the reading of the BCR issue at the transfer station