

Title:	Export not finished	ID:	
		0104	
Date in:	Response:	Model:	Author:
2003-03-29	2003-03-29	-	CMa

Q:

I have a problem here with one of the robots.

Here's the sequence of commands that I've issued.

st 1900

OK

st 1801

OK

rd 1915

0

rd 1915

1

wr dm0 4

OK

wr dm5 20

OK

st 1905

OK

and then the robot gets the plate from the hotel and then it stops.

i issued

rd 1814

0

but

rd dm200

09728

rd 1915

0

The mover has no mechanical problem. It was verified by your friend that was here.

Why is there no error reported? flag in 1814 should be set to 1. Can you look into this problem? So far I've seen this problem in two of the robots and I haven't checked one of them yet.

Also, one of the robot is moving very slowly. It's making loud sounds as the robot goes up and down. Do you have any ideas as to what might cause this problem?

A:

Urs mentioned something that you had difficulties one of your computers. Any news on this?

Check if input 102 or 103 is on. The system could assume there's a active transfer station or try to continue the export with ST 1901.