

| | | | |
|-----------------|--|---------------|----------------|
| Title: | How to detect the glass door switch | ID: | |
| | | 0133 | |
| Date in: | Response: | Model: | Author: |
| 2003-09-19 | 2003-09-19 | - | CMa |

Q:

Can you please tell me how to read the Glass door sensor?
Also, is there any way to query the device to find out if it has a user gate or swap station?
If possible, I'd like an exhaustive list of the possible commands that I can use to acquire as much information from the system as possible.

A:

User Door : RD 1811 (in ActiveX we poll 1814, 1811 and 1915)
Swap Stn: RD 3210 ('1' for active = motorized x-fer Stn.)

There is no such list. But feel free to ask. I'm happy to forward you the information.