

| | | | |
|-----------------|----------------------------|---------------|----------------|
| Title: | Unit not responding | ID: | |
| | | 0137 | |
| Date in: | Response: | Model: | Author: |
| 2005-03-08 | 2005-03-08 | - | WOp |

Unit not responding

1. check connection from PC to the unit (correct connectors and cables)
2. check the communication settings
 - 1) 9600 Baud
 - 2) 8 data bits
 - 3) 1 stop bit
 - 4) Parity even
3. make sure unit is powered ON