

Title:	Shovel potentiometer adjustment no reaction			ID:
				0192
Date in:	Response:	Model:	Author:	
2004-08-22	2004-08-22	HS Carousel	CMa	

Q:

There is no reaction on the potentiometer adjustments for the shovel. This is a HS-system.

A:

Just talked with Yuri. To me it looks like T48 has not been updated. T48 is the amount of time the shovel runs at its full speed. You can read this timer by "RD T48". The last number holds the time. By "WS T48 xx" you can permanently change this timer. You may need to shorten this time. E.g. IF "RD T48" gives "0.00000,00045" you may change it by "WS T48 35".

Note that the potentiometer settings will only be valid after T48 has elapsed.