



Title:	Carrousel Connector Layout			ID:
				0021
Date in:	Response:	Model:	Author:	
2004-06-29	2004-06-29	STX200	CMa	

Q:

We have a problem with the carrousel positioning. Can you send me a wiring diagram?

A:

Please find attached data..

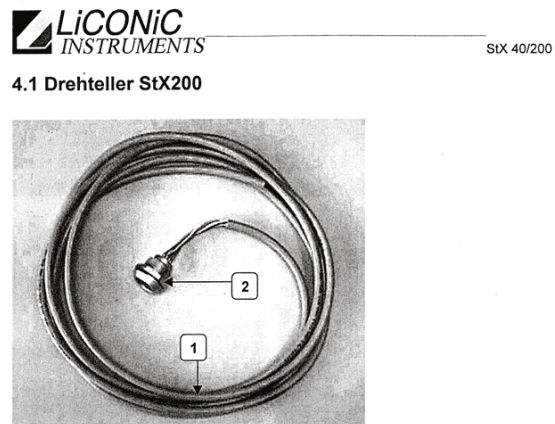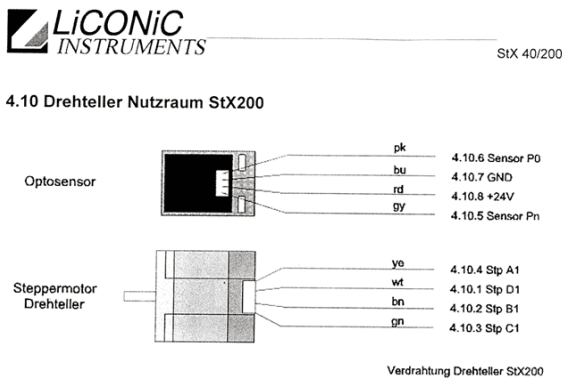

Kabeldefinition

Pos	Name	Stk/Dim	Bezeichnung	Lieferant	Art-Nr.
1		2500mm	Kabel 8x0.5		
2	4.1		Buchse Lemo 8polig ERA.2E		

Kabelbelegung

Stecker/Pin	Farbe/Nr.	Kabel Pos	Bezeichnung	Stecker/Pin
4.1.1	wt	1	Stepper D1	1.1.10
4.1.2	bn	1	Stepper B1	1.1.9
4.1.3	gn	1	Stepper C1	1.1.8
4.1.4	ye	1	Stepper A1	1.1.7
4.1.5	gy	1	Sensor Pn	1.1.2
4.1.6	pk	1	Sensor P0	1.1.1
4.1.7	bu	1	Sensor GND	1.1.5
4.1.8	rd	1	Sensor +24V	1.1.3