

<b>Title:</b>	<b>Swap Station 1912 command does not work</b>			<b>ID:</b>
				0280
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>	
2007-05-26	2007-05-26	-	CMa	

## Q:

We replaced the standard TurnStation against a SwapStation. The 1912 command does not work.

## A:

The SwapStation works in SlaveMode (the user defines the station position). The SlaveMode must be activated in the FW. You need to install a new FW with the SlaveMode (3105) activated.