

Title:	Activate Shovel Service Mode	ID:	
		0003	
Date in:	Response:	Model:	Author:
2006-03-31	2006-03-31	-	CMa

Q:

My system stopped at the stacker in an Unload. How can I activate the shovel to check if the the shovel is positioned correctly?

A:

Here is the sequence to move the shovel:

FW Version <=6x

```
WR DM123 63450 ; Activate service mode
ST 1701        ; shovel out
RS 1701        ; shovel in
WR DM123 0     ; Leave service mode
```

FW Version 7ff

```
WR DM123 63450 ; Activate service mode
ST 1401        ; shovel out
RS 1401        ; shovel in
WR DM123 0     ; Leave service mode
```