

Title:	Black Knob Functionality for Carrousel	ID:	
		0360	
Date in:	Response:	Model:	Author:
2010-06-18	2010-06-18		CMa

Q:

When the multi-function switch on the front panel is depressed to reset an alarm, it also causes the carousel to index several degrees. If we then resume a test which was already in progress, the device does not recognize that the stacker had been displaced from its original position, and the handler goes to pick the wrong plate. This behaviour is really undesirable.

Questions:

1. What are the intended functions of this switch/
2. What can be done to prevent the behavior I described above?

A

When system is initialized and in idle (1903) the switch does

Short press -> move one position

Long press -> move 1/3 rev.

This is to look at the unit's content without need to open.