

Title:	Cooler LEDs			ID:
				0363
Date in:	Response:	Model:	Author:	
2016-12-24	2016-12-24		CMA	

Cooler LED's

LED	Perm.	Blinking Slow	Blinking Fast	Blinking Intermediate
red	No Power-Voltage	General Error	Warning Rest	
yellow	1 st Stage ON	1 st Stage Enabled	1 st Stg. OverTemp / OFF	1stOverPress
green	2 nd Stage ON	2 nd Stage Enabled	2 nd Stg. OverTemp / OFF	2ndOverPress Stepper Valve Moving