

Title:	Stepper Valve Positioning			ID:
				0393
Date in:	Response:	Model:	Author:	
2014-04-17	2014-04-17		CMA	

Q:

Stepper Valve position was set to „0“

A:

The stepper valve should not be on “0”. This would mean that it is completely closed. If suspicious you may re-initialize the stepper valve by switching the cooler off and restart. You can hear the initialization process if pressing your ear against the housing. The initialization is time based, since the valve has no end-switch. The homing movement is 30s fixed. When reaching the home position you can hear a clicking. Next you'll hear a short stop and then the actual positioning.