


Title:	PLC System Reset			ID:
				130
19.09.24	Response:	Model:	Author:	
		-	OKr	

PLC Factory Reset


Issue: PLC blocked, with display red



- Step 1:
- Remove power
 - Press and hold up and down key simultaneously
 - While holding the keys down, power on the unit and wait for the display to flash green

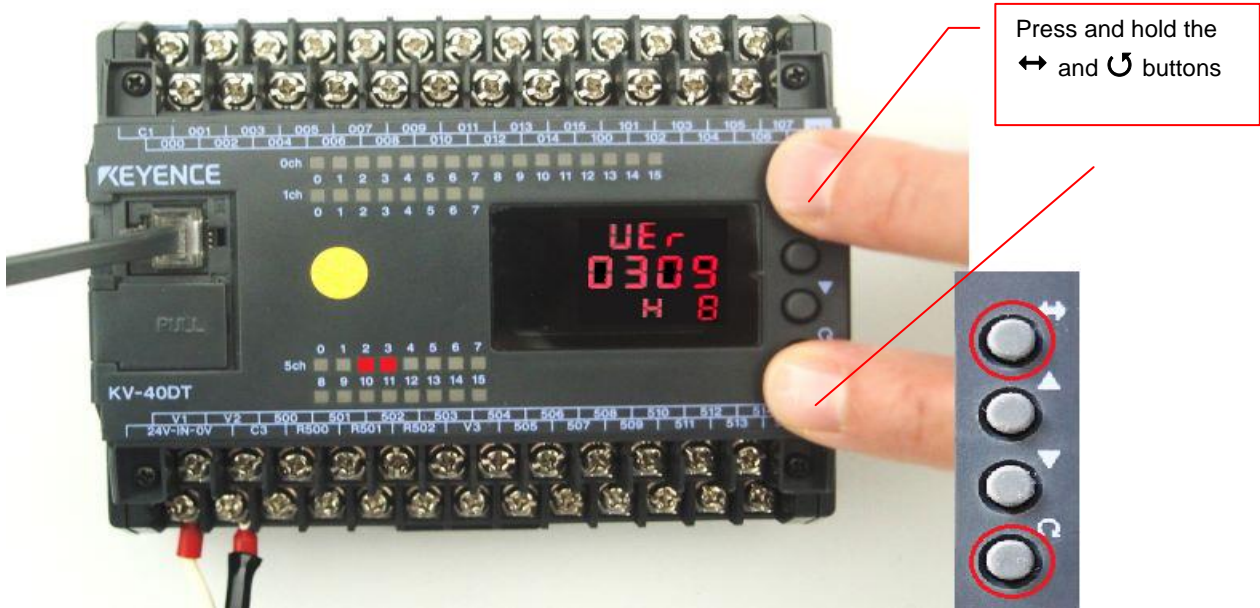


- Step 2: - Press and hold ← and ↻ simultaneously
 - A green flashing test pattern will appear on the display

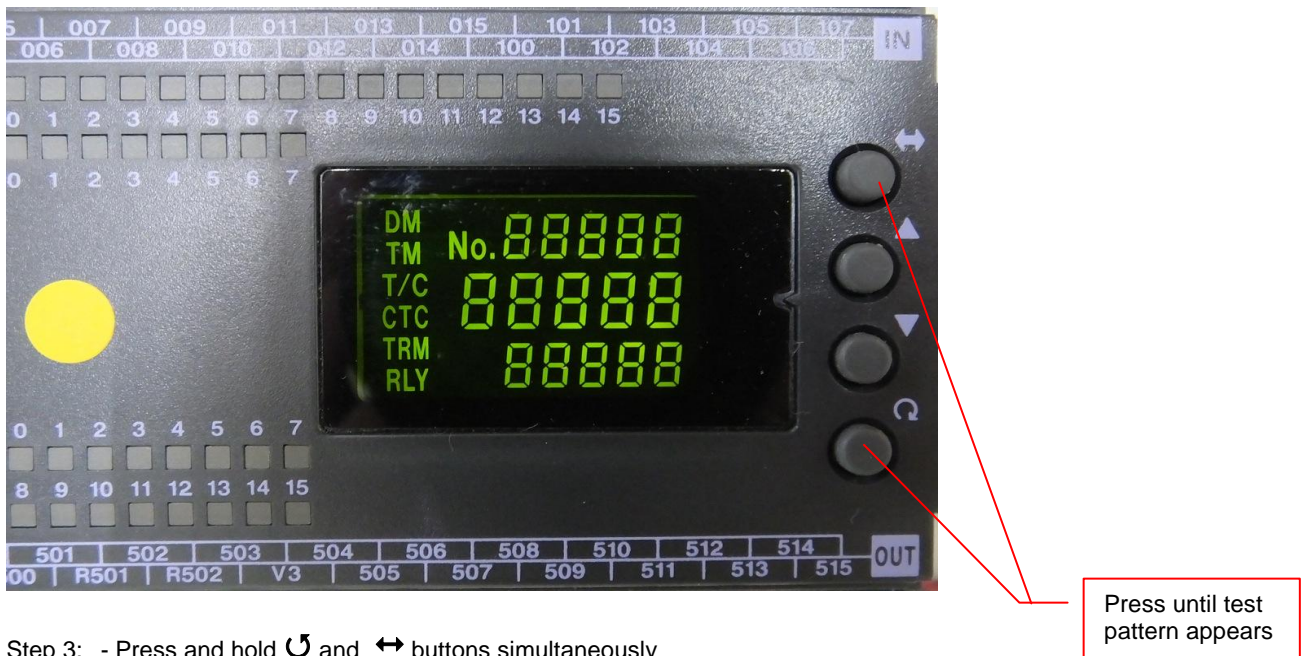


- Step 3: - Press and hold ↻ and ← buttons simultaneously
 - Wait for flashing to stop

- Step 4 - Push the ↻ and ↔ buttons simultaneously again
 - Wait for the display to flicker between green and red

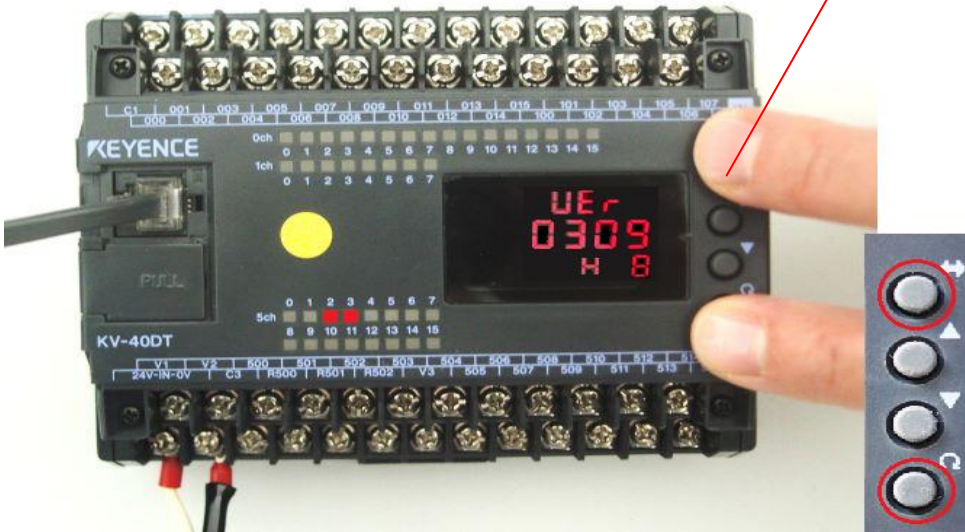

after pushing twice you will get this result on the display

- Step 5 - Push the ↻ and ↔ buttons simultaneously again
 - Wait for above message to appear on the display

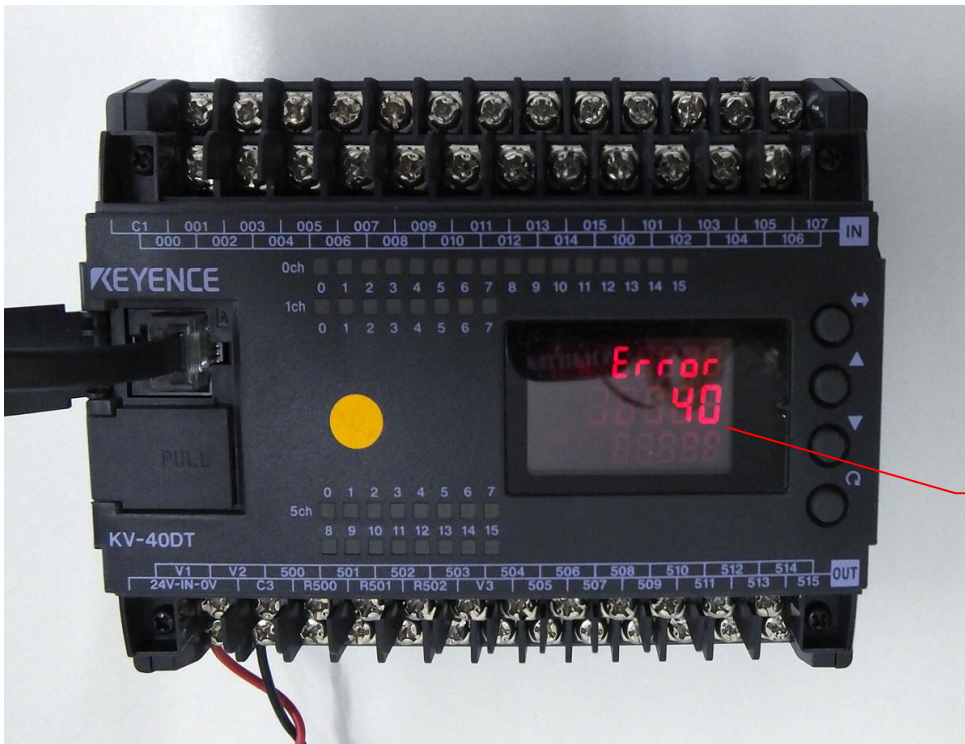

after you turn power back on, this should appear

- Step 6: - Power off PLC and wait 10 seconds
 - Turn the power back on
 - The text "Error 40" should now appear on the display
 - The PLC is now ready for re-programming