


Title:	Flush Gate Motors			ID:
				148
30.11.16	Response:	Model:	Author:	
		-	CMA	

Flush-Gate Motors

There three versions of Flush-Gate motor-gearbox combination: V2.1, V2.2, V2.3. The gearboxes have some incompatibilities

- ! Diameter of axis V2.1: 8mm, V2.2 and V2.3: 10mm
- ! Attachment holes and outer dimensions V2.1 differ from V2.2 and V2.3

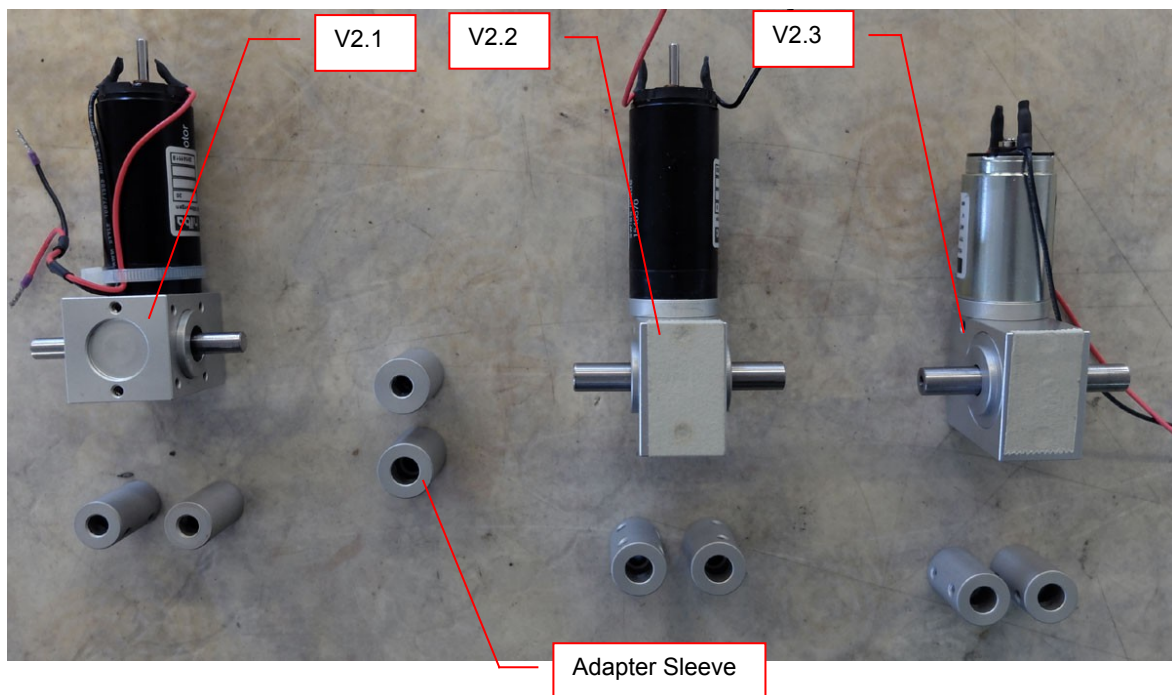

Version 2.2 and V2.3 are interchangeable. To replace a version 2.1 motor-gearbox two Adapter Sleeves are required.