

Title:	<b>Shaker Cassette Optimization</b>			ID:
				0015
Date in:	Response:	Model:	Author:	
2007-05-09	2007-05-09	-	C <i>Ma</i>	

## Fine Tune Shaker Cassettes

On high shaking speeds or amplitudes plates may slide out of the cassettes. The retaining properties may be optimized by:

- Retaining Lip Height to be 1 .. ~1.5 mm
- Retaining Lip on left and right Side Wall to be identical
- Minimize play of plate by optimizing Center Lip
- Make undercut at the end of Plate Rail



