


Title:	Condensation on Glass Door			ID:
				0090
Date in:	Response:	Model:	Author:	
2010-05-05	2010-05-05	-	CMa	

Condensation on Inner Glass Door

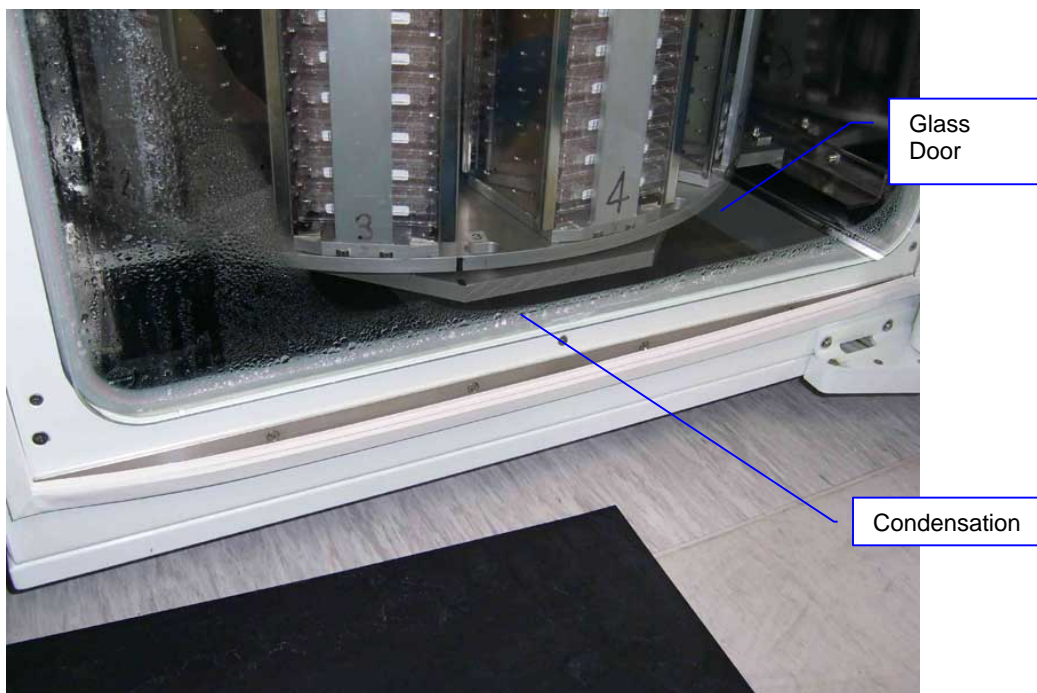

I assume that the condensation is on the inner side of the glass door and the unit is set to high humidity. Basically there are two reasons for excessive condensation known.

1) Temperature between outer door and glass door too low: Are the door heater working? Do they open the outer door often or leave it open?

2) RH control: Pls. check the RH settings (RD DM893). Could be that the RH has to be re-calibrated because it is too high.